twinnno

RM405 手持拉曼分析儀
Handheld Raman Analyzer

使用說明書 **User Manual**

Ver. 1.0

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Table of Contents

1 Introduction	23
1.1 Description of Contents	23
1.2 Exterior Description	24
1.3 Product Features	25
1.4 Instructions for Product Use	26
2 Basic Uses	27
2.1 Power On/Power Off	27
2.2 Micro SD Card Installation Method	28
2.3 Measurement Method	29
2.4 System Reset	30
2.5 Battery Charging Mode	31
2.6 USB Mode.....	32
3 System Operation	33
3.1 MATCH Measurement Description	34
3.2 COMP Measurement Description	36
3.3 Setup Operating Instructions	37
3.4 BROWSER Operating Instructions	38
3.5 SYSTEM Operating Instructions	39
4 Specification	40
5 Limited Product Warranty.....	41

1 Introduction

1.1 Description of Contents

The shipment should contain the following objects. If anything is missing, please contact the distributor.


Suitcase

RM405

**Transmission
Line**

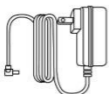Charger

Extraction Bottle
*6



Warranty Card

Manual

1.2 Exterior Description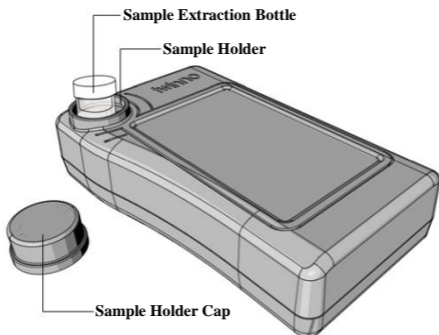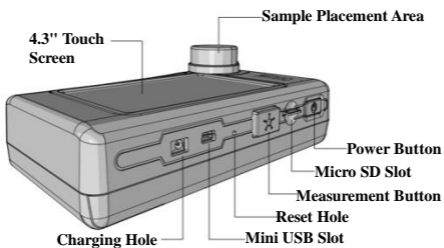

1.3 Product Features

1. Built-in 785nm laser.
2. Internal 405nm .
3. Wide range Raman shift ($400\text{cm}^{-1} \sim 1.4000\text{cm}^{-1}$).
4. Support SERS measurement.
5. Built-in Raman spectra database, also can 1. create database by user.
6. Support data IO and compare functions.
7. Nondestructive analysis with rapid measurement capability.
8. The measurement procedure is simple, without complex pretreatment.
9. Simply and easily operation.
10. System measurement design for stand-alone without computer.

1.4 Instructions for Product Use

1. RM405 is high precision optical measurement equipment. Any collision, vibration or fall can affect measurement accuracy.
2. To assure measurement accuracy, we suggest send the unit back to us for calibration once per year.
3. RM405 uses lithium battery. Please do not store or operate it under high temperature condition to prevent the battery from combustion or explosion.
4. During charging, if overheating, smoking or odor occurs, please stop charging immediately.
5. Please do not remove or alter the battery.
6. During charging, do not place the equipment in hot environment, do not use cloth or others that affect heat dissipation to cover or wrap it. Otherwise, it may cause equipment fault due to poor heat dissipation.
7. Please do not use organic solvent to clean the equipment. Otherwise, it may damage the product exterior or touch screen.
8. If the product cannot be used or operated normally, please do not disassemble and repair it literally.
9. Please do not use hand rope or similar thing to hang or swing the equipment to prevent damage.
10. Product function change without notice, please check at the official website to update.


2 Basic Uses

2.1 Power On/Power Off

Power On: Short press the power button, green light will go on.

Power Off: Long press the power button for 3 seconds, green light will go off.


Diagram for reference: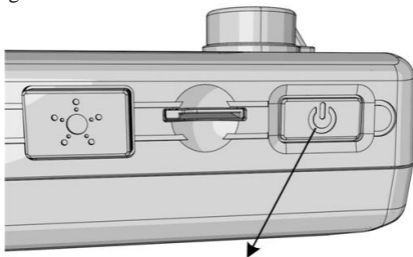


**Power on will have the green light on.
Power off will have the green light off.**

2.2 Micro SD Card Installation Method

Install Micro SD card: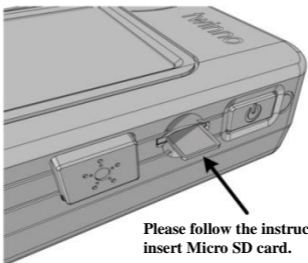

Take out Micro SD card:

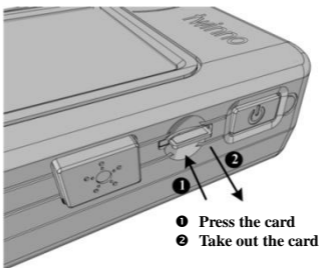2.3 Measurement Method

Process flow:

1. Put liquid sample into the extraction bottle.
2. Tighten the extraction bottle cap.
3. Put extraction bottle into the sample holder.
4. Put on the sample holder cap and tighten it.
5. Press the measurement button and start measurement.
6. Speaker sounds and the upper right corner Standby indicates measurement completed.


Precautions:

1. During measurement, we suggest to put on the cap to reduce the impact from environment light.
2. Prior to using extraction bottle, please first clean dust inside.
3. Keep cleanness on extraction bottle surface. Wipe clean if fingerprint is found.
4. The liquid level shall be higher than a half the extraction bottle.
5. The liquid level in the extraction bottle shall be higher than the minimum required level.
6. During measurement, please place it horizontally, and prevent shaking and tilting the liquid level from affecting the measurement accuracy.

2.4 System Reset

When the system is down, please press reset button to reset the system.

Diagram for reference:

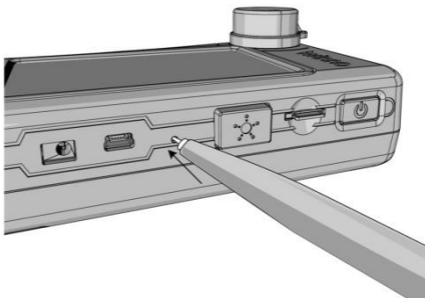Suggest to use pencil tip pressing to reset the system

2.5 Battery Charging Mode


1. Connect charger to RM405 charging hole to start charging.
2. For charging when power is off, power button will show red light. After charging is completed, red light will go off.
3. For charging when power is on, the upper right corner of screen will show a battery logo and the power button will show red light. After charging is completed, red light will go off.

Precautions: RM405 does not provide USB charging function.

Diagram for reference:

During charging the power button will show red light. When charging is completed, the red light will go off.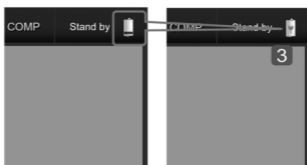

During charging, battery power level symbol will be changed to charging status symbol

2.6 USB Mode

1. RM405 supports USB. Users can select storage device mode or PC connection mode.
2. Storage device mode (MSC): access micro SD.
3. PC connection mode (SPD): use Opticker program to operate RM405 measurement from PC.

Diagram for reference:


USB has SPD and MSC modes for selection

Screen after connecting USB transmission cable


3 System Operation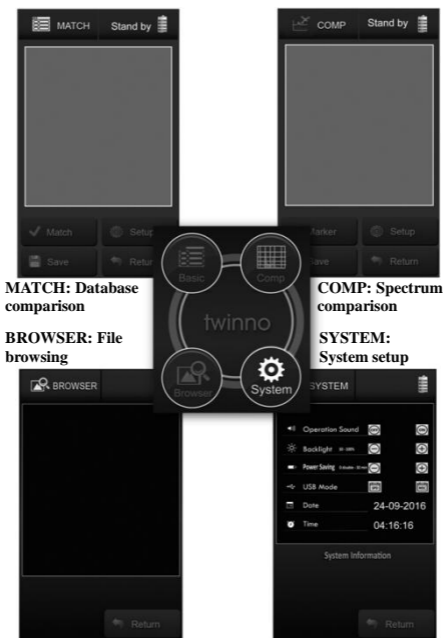


3.1 MATCH Measurement Description

Description: compare Raman signal to database, identify material composition.


MATCH interface: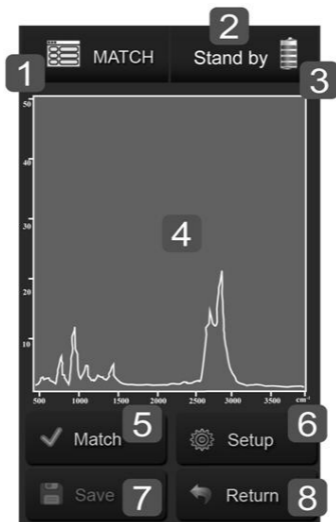

1. System mode.
2. System status.
3. Battery power level.
4. Measurement result (Raman spectrum).
5. Comparison function -- after measurement press the button to indicate whether the sample is a substance in the database. If the database does not have the substance, it is not identified, as in the following figure.
6. Parameter setup.
7. Store measurement data.
8. Return to previous page.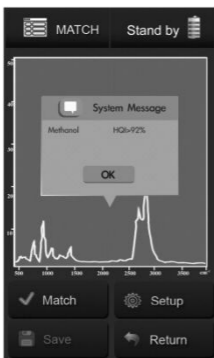

Database comparison result.

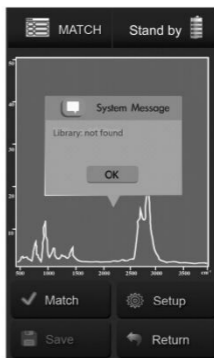
Database does not found. It cannot be identified.

3.2 COMP Measurement Description

Description: Simultaneously show several measurement results on the same coordinate axis for user comparison.

COMP interface: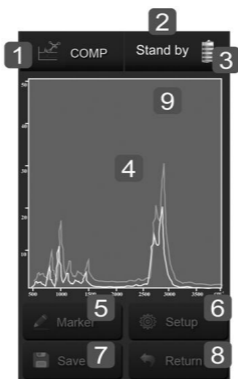

1. System mode.
2. System status.
3. Battery power level.
4. Measurement result (Raman spectrum).
5. Record measurement spectrum.
6. Parameter setup.
7. Store measurement data.
8. Return to previous page.
9. Spectrum color indication -- press here to clear the spectrum on screen.

3.3 Setup Operating Instructions

Description: Users can set up measurement parameters by themselves.

Diagram for reference:


1. Measurement method.
2. Exposure time (10 ~ 50000ms).
3. Laser power (10 ~ 100%).
4. Laser wavelength calibration.
5. Smooth wave filtering parameter.
6. Average times.
7. HQI parameter.
8. Return to previous page.

3.4 BROWSER Operating Instructions

Description: Users can view the measurement data (File) on micro SD card and database (Library).

BROWSER interface: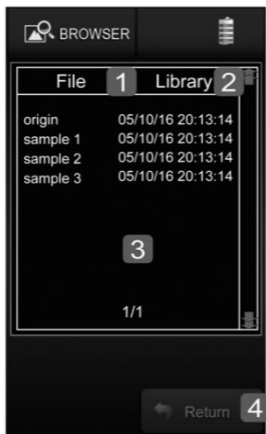

1. Measurement data.
2. Database.
3. File name information.
4. Return to previous page.

3.5 SYSTEM Operating Instructions

Description: Users can set up system parameters by themselves

SYSTEM interface:

1. Sound setup (ON, OFF).
2. Backlight setup (30% ~ 100%).
3. Automatic power off time setup (0 ~ 30 minutes).
4. USB mode setup (MSC, SPD).
5. Date setup.
6. Time setup.
7. System information (hardware version, firmware version, optical drive serial number, Device license).
8. Return to previous page.

4. Specification

Sensor	Raman shift range	400cm ⁻¹ ~ 3000cm ⁻¹
	Exposure time	10ms ~ 50s
	Resolution	~25cm ⁻¹
Laser	Excited wavelength	785nm ±5nm
Measure-ment	Mode	1. MATCH: Substances Detection 2. COMP: Signal Comparison 3. BROWSER: Data Browser 4. SYSTEM: System Configuration
Other	Input/output interface	micro SD (SDHC maximum 32GB), USB2.0
	Data format	.jmr
	Battery	2600 mAh rechargeable lithium battery ×1
	Battery usage time	Maximum 4 hours (fully charged)
	Charger	5V/2A
	Screen	4.3" LCD (480 × 800 pixel), including touch control function
	RM405 dimensions	155mm × 95mm × 60mm (L × W × H)
	Extraction bottle dimensions	15mm × 50mm (Dia. × H)
	Weight	500g
	Operating temperature	15 ~ 30°C
	Storage temperature	0 ~ 40°C
Language	English	

5. Limited Product Warranty

strider tech guarantees that RM405 products will provide consumers with free repair or replacement for the products of the same grade during the warranty period, such as defects or malfunctions in materials or functions. This warranty does not cover any damage resulting from improper installation, accident, abuse, misuse, natural disaster, inadequate or excessive power supply, abnormal mechanical or environmental conditions, and any unauthorized removal, repair or modification of the product.

The warranty does not apply to the following situations:

1. Users do not follow the instructions to use the product.
2. Users do not use the product for its designed use. If any original identification information on the product has been altered, erased or removed, or if any product has not been properly handled or packed, sold as a second-hand item, or resold in violation of local export laws, the limited product warranties will be void.

This limited product warranty covers only the repair, replacement, or refund of the defective product as described above. Our company shall not be liable for any damages or loss resulting from the loss, damage, or error of the contents or data, or for determining the source of the system's problem, or the demolition, repair or installation of the product, and nor is it covered by the product warranty. This warranty does not cover third-party software, connected devices, or stored data. Therefore, we shall assume no responsibility for any loss or damage caused by the third party's software, connected equipment or stored material.

5.1 Product within Warranty Period

1. RM405 has 2 years of original manufacturer warranty.
2. Consumables for the product are not covered by the warranty.
3. New product replacement:

Consumers within 7 days after the purchase of a product that has functional defects or failure, can apply for a replacement.

Note: The returned product must be in the original packaging with all parts included.

4. Maintenance services:

For product that has functional defects or failure within 7 days of purchase, the original manufacturer provides free repair.

Note: After receiving the returned product, the original manufacturer will conduct basic check. If the check result finds it is not the product fault and failure, the warranty will not apply.

5. How to send for repair:

Users return the product to the agent or directly to the original manufacturer.

5.2 Product is Outside the Warranty Period


After the warranty period has passed for more than 2 years, Users can still return the product of fault or malfunction to the original manufacturer for repair. Our company will charge the users for repair. If the model is discontinued, we have the right to replace it with an equivalent new product.

Opticker Download And Install

1. Connect to **twinno** official website :

www.twinno.com

2. 

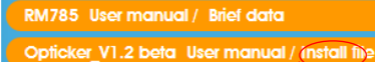3. 

Click Install file

4. Download file and install Opticker_Install.exe.
5. After install, Opticker.exe path appear on desktop.
6. Execute Opticker.exe start up the program.

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